

Niton XL3 Analyzer Quick Start Guide

IMPORTANT SAFETY INFORMATION:
We recommend that you read the entire user's manual before using your analyzer

1




Connect the Charger To the Wall Outlet



Orient the Battery Pack Properly

MAKE SURE THAT THE KEY IS IN FRONT AND THE THIN SIDE IS DOWN

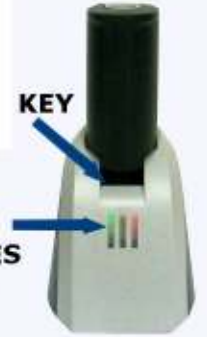

GREEN LIGHT INDICATES FULLY CHARGED

Place Battery in Charger

2



PUSH BATTERY PACK INTO HANDLE, THIN SIDE UP, THEN SECURE LATCH

Insert Battery



Analyzer Controls

ON/OFF/ESCAPE BUTTON

4-WAY TOUCH PAD

CLEAR/ENTER BUTTON

INDICATOR LIGHTS

OPEN COVER TO ACCESS PORTS


PRESS AND HOLD ON/OFF/ESCAPE BUTTON UNTIL TOUCH SCREEN POWERS ON



Power Up

3

TOUCH SCREEN

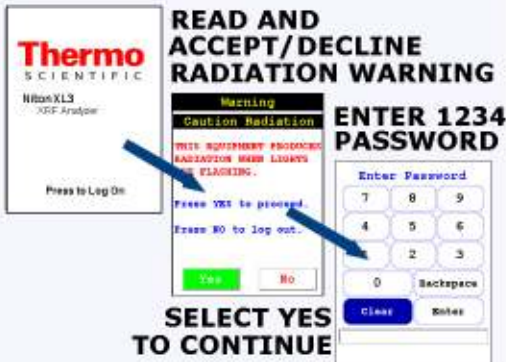

Startup Screens



Perform System Check




Main Menu
PART NUMBER 410-008

Niton XL3 Analyzer Quick Start Guide

IMPORTANT SAFETY INFORMATION:
We recommend that you read the entire user's manual before using your analyzer


4

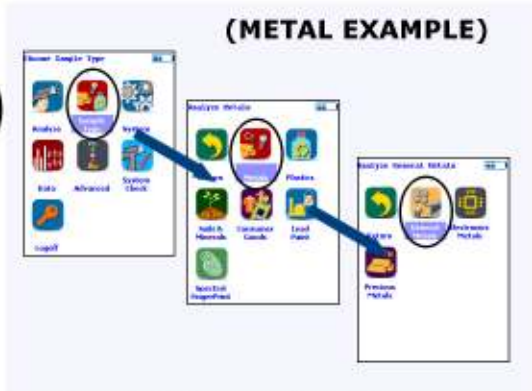

Select Sample Type For Analysis




From Ready To Test Screen, Analyze Alloy Check Sample Provided

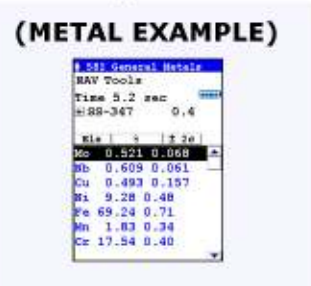

Stop Reading After 10 Secs

REFER TO THE USER'S GUIDE FOR SPECIFIC STANDARD OPERATING PROCEDURE


5

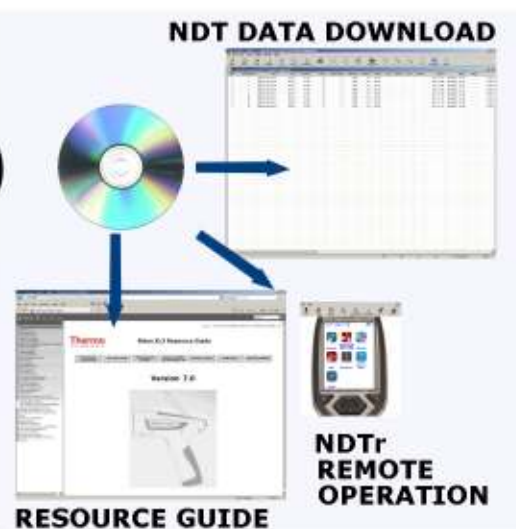

Install PC Software Resource Guide/ User's Guide



Start PC Software

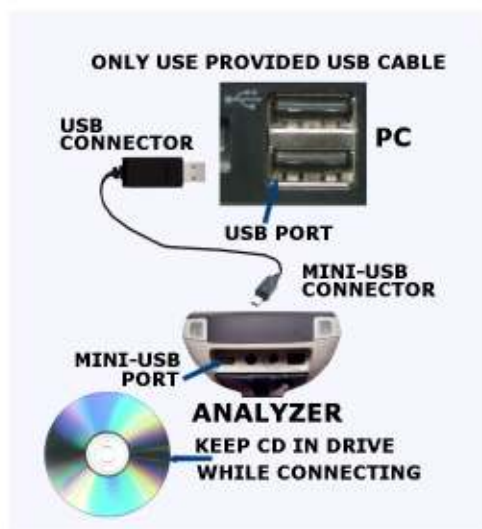


Install USB Driver Provided On Disk

6



Connect to PC

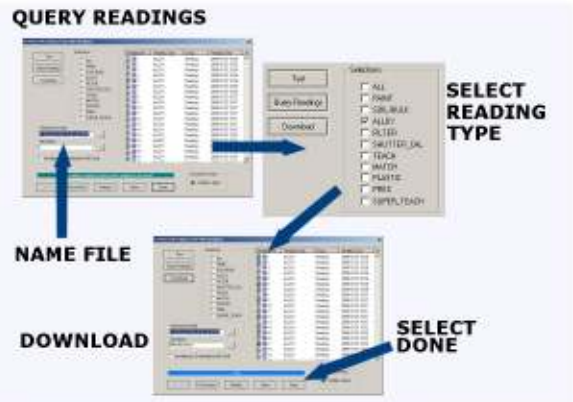

Download Data to PC



Data Report